

Cytek® cFluor® Immunoprofiling Kit

PRODUCT DETAILS																																																															
Catalog number:	R7-40000 (25 Tests)																																																														
Category:	Immunophenotyping																																																														
Format:	cFluor® conjugated antibodies in individual tubes																																																														
	<table> <tr> <th>Target</th><th>Clone</th><th>Fluorochrome</th><th>Item number</th></tr> <tr><td>CD3</td><td>SK7</td><td>cFluor® V420</td><td>R7-20054</td></tr> <tr><td>CD4</td><td>SK3</td><td>cFluor® R780¹</td><td>R7-20084</td></tr> <tr><td>CD8</td><td>SK1</td><td>cFluor® B515²</td><td>R7-20036</td></tr> <tr><td>CD14</td><td>M5E2</td><td>cFluor® V450</td><td>R7-20004</td></tr> <tr><td>CD16</td><td>3G8</td><td>cFluor® R668²</td><td>R7-20070</td></tr> <tr><td>CD19</td><td>H1B19</td><td>cFluor® BYG710¹</td><td>R7-20010</td></tr> <tr><td>CD25</td><td>BC96</td><td>cFluor® BYG781¹</td><td>R7-20080</td></tr> <tr><td>CD27</td><td>QA17A18</td><td>cFluor® R840^{1,2}</td><td>R7-20082</td></tr> <tr><td>CD45</td><td>HI30</td><td>cFluor® V547²</td><td>R7-20012</td></tr> <tr><td>CD45RA</td><td>HI100</td><td>cFluor® B690¹</td><td>R7-20086</td></tr> <tr><td>CD56</td><td>5.1H11</td><td>cFluor® R720²</td><td>R7-20090</td></tr> <tr><td>CD127</td><td>A019D5</td><td>cFluor® R659</td><td>R7-20078</td></tr> <tr><td>CD197 (CCR7)</td><td>G043H7</td><td>cFluor® BYG575</td><td>R7-20076</td></tr> <tr><td>IgD</td><td>IA6-2</td><td>cFluor® BYG667¹</td><td>R7-20088</td></tr> </table>	Target	Clone	Fluorochrome	Item number	CD3	SK7	cFluor® V420	R7-20054	CD4	SK3	cFluor® R780 ¹	R7-20084	CD8	SK1	cFluor® B515 ²	R7-20036	CD14	M5E2	cFluor® V450	R7-20004	CD16	3G8	cFluor® R668 ²	R7-20070	CD19	H1B19	cFluor® BYG710 ¹	R7-20010	CD25	BC96	cFluor® BYG781 ¹	R7-20080	CD27	QA17A18	cFluor® R840 ^{1,2}	R7-20082	CD45	HI30	cFluor® V547 ²	R7-20012	CD45RA	HI100	cFluor® B690 ¹	R7-20086	CD56	5.1H11	cFluor® R720 ²	R7-20090	CD127	A019D5	cFluor® R659	R7-20078	CD197 (CCR7)	G043H7	cFluor® BYG575	R7-20076	IgD	IA6-2	cFluor® BYG667 ¹	R7-20088		
Target	Clone	Fluorochrome	Item number																																																												
CD3	SK7	cFluor® V420	R7-20054																																																												
CD4	SK3	cFluor® R780 ¹	R7-20084																																																												
CD8	SK1	cFluor® B515 ²	R7-20036																																																												
CD14	M5E2	cFluor® V450	R7-20004																																																												
CD16	3G8	cFluor® R668 ²	R7-20070																																																												
CD19	H1B19	cFluor® BYG710 ¹	R7-20010																																																												
CD25	BC96	cFluor® BYG781 ¹	R7-20080																																																												
CD27	QA17A18	cFluor® R840 ^{1,2}	R7-20082																																																												
CD45	HI30	cFluor® V547 ²	R7-20012																																																												
CD45RA	HI100	cFluor® B690 ¹	R7-20086																																																												
CD56	5.1H11	cFluor® R720 ²	R7-20090																																																												
CD127	A019D5	cFluor® R659	R7-20078																																																												
CD197 (CCR7)	G043H7	cFluor® BYG575	R7-20076																																																												
IgD	IA6-2	cFluor® BYG667 ¹	R7-20088																																																												
Test Dilution:	5µl per test																																																														
Application:	Flow Cytometry																																																														
Formulation:	Phosphate-buffered saline, pH 7.2, containing 0.09% sodium azide and 0.2% BSA (BSA Country of Origin)																																																														
Storage:	2-8°C and protected from light Do not freeze																																																														

PRODUCT DESCRIPTION

The Cytek® cFluor® Immunoprofiling Kit allows for the identification of helper T cells, cytotoxic T cells, B cells, NK cells, and monocytes in human peripheral blood mononuclear cells and in whole blood. The reagents in this kit help to distinguish different subsets of T, B, and NK cells; including regulatory T cells, naïve T cells, activated T cells, memory T cells, effector T cells, naïve B cell, memory B cells, nonclassical, and classical monocytes.

CD3 is expressed on all mature T cells, NKT cells, and some thymocytes. CD3, a part of the CD3/T cell receptor complex, plays a role in antigen recognition, signal transduction, and T cell activation.

CD4, is expressed on most thymocytes, a major subset of T cells, and on monocytes/macrophages. Functionally, CD4 is associated with thymic differentiation, in conjunction with MHC class II molecules in antigen recognition and with signal transduction.

R8-50000 Rev. C, Effective Date: 03/17/2025

Cytek Biosciences Inc. | 47215 Lakeview Boulevard, Fremont, CA 94538

Phone: +1 (510) 657-0102 | Fax: +1 (510) 657-0151 | cytekbio@cytekbio.com | www.cytekbio.com

CD8 is found on thymocytes, on a subset of T cells, and on NK cells. This molecule acts as a co-receptor in MHC class I molecules in antigen recognition and has a role in T cell activation and in thymic differentiation.

CD14 is also known as a high affinity LPS receptor and is highly expressed on monocytes and macrophages. It is also expressed on granulocytes, but at a lower level. In addition, CD14 is found on interfollicular dendritic cells, reticular dendritic cells, and Langerhans cells.

CD16 is expressed on NK cells, monocytes and macrophages in the form of CD16a. Another form of CD16, CD16b is expressed on neutrophils. CD16 engagement of IgG leads to NK cell activation, antibody-dependent cell-mediated cytotoxicity (ADCC) and phagocytosis.

CD19 is expressed in the B cell lineage from pro-B to blastoid B cells. However, it is absent on plasma cells. It is also expressed on follicular dendritic cells. CD19 is involved in B cell development, activation, and differentiation.

CD25 is a low affinity IL-2 receptor that is expressed on progenitor lymphocytes, activated T cells, B cells, and monocytes, as well as on regulatory T cells. When CD25 is in association with CD122, IL-2 receptor beta chain, and with CD132, the common gamma chain, the resulting complex is the high-affinity IL-2 receptor. IL-2 receptor signaling is key lymphocyte proliferation and survival.

CD27 is a lymphocyte specific member of the TNF receptor superfamily. CD27 is expressed on a subset of thymocytes, mature T cells. The expression further upregulated in T cell activation. Subpopulations of B cells and NK cells also express CD27. In T cell-B cell interaction, CD27 binds to CD70 to provide costimulation to result in T cell activation and B cell differentiation and proliferation.

CD45 is expressed on all hematopoietic cells, except erythrocytes and platelets. CD45 is a signaling molecule that is involved in cellular proliferation, differentiation and in regulation of immune cell functions.

CD45RA is an isoform of CD45 due to alternate splicing. CD45RA is expressed on naïve or resting CD4+ and CD8+ T cells, as well as on B cells and monocytes.

CD56, within the hematopoietic system is expressed on NK cells and NKT cells, a subset of T cells. In the nervous system, CD56 is expressed by neurons and plays a role in the homotypic adhesion of neural cells.

CD127, when it heterodimerizes with the common gamma chain (CD132), forms the IL-7 receptor. This receptor is involved in the development of B cells in the bone marrow, and the control of T cell proliferation in the periphery. CD127 expression is downregulated in regulatory T cells.

CD197 (CCR7) is a chemokine receptor for chemokines CCL19 and CCL21. CCR7 and its ligands are involved in trafficking and migration of immune cells into immunological milieus. For example, B cell migration and organization into secondary lymphoid tissues, and migration of memory T cells into inflamed tissue. CCR7 is expressed on T cells and the detection of CCR7 and, in combination with CD45RA, can be used to distinguish naïve T cells from central and effector memory T cells.

IgD is a molecule that is expressed on naïve B cells. IgD is expressed on B cells upon exiting the bone marrow to the periphery.

REFERENCES

Clevers H et al. Ann Rev Immunol. 1988 **6**, 629 (1988)
 Evans RL et al. Immunol. **78**, 544 (1981)
 Cole DK et al. Cell Mol Immunol. **1**, 81 (2004)
 Ferrero E et al. J. Immunol. **145**, 331 (1990)
 Wirthmueller U et al. J Exp Med. **175**, (1992)
 Dennis O et al. J Immunol. **170**, 73 (2003)
 Shevach E. Nat Rev Immunol. **2**, 389 (2002)
 Vossen M et al. J Immunol. **180**, 3739 (2008)
 Fritsch R et al. J Immunol. **175**, 6489 (2005)
 Loken MR et al. Cytometry. **11**, 453 (1990)
 Kumar, Connors & Farber. Immunity. **48**, 202 (2018)
 Poli A et al. Immunology. **126**, 458 (2009)
 Carrette, F & Surh, C. Semin Immunol. **24**, 209 (2012)
 Mahnke Y et al. Eur J Immunol. **43**, 2797 (2013)
 Goodwin R et al. Cell. **73**, 447 (1993)

RECOMMENDED USAGE

Human peripheral mononuclear cells (PBMC) or whole blood has been tested to validate the performance of this kit. Please refer to the product web page for the staining protocol, fluorochrome tag list, experiment template, and data acquisition protocol.

Please briefly centrifuge the reagent vial before use.

Use appropriate personal protective equipment per the product safety data sheet when using this product.

¹cFluor® BYG667, cFluor® BYG710 and cFluor® BYG781 are tandem dyes made with R-PE. cFluor® B690 is a tandem dye made with PerCP. cFluor® R780 and cFluor® R840 are tandem dyes made with APC. Caution – Tandem dyes may show changes in their emission spectra with prolonged exposure to light or fixatives.

²cFluor® B515, cFluor® R668, cFluor® R720, cFluor® R840, and cFluor® V547 are equivalent to CF®488A, CF®647, CF®700, APC-CF®790T, and CF®405L respectively, manufactured and provided by Biotium, Inc. under an Agreement between Biotium and Cytek (LICENSEE). The manufacture, use, sale, offer for sale, or import of the product is covered by one or more of the patents or pending applications owned or licensed by Biotium. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. No right under any other patent claim, no right to perform any patented method, and no right to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is conveyed expressly, by implication, or by estoppel.

"Cytek" and "cFluor" are trademarks of Cytek Biosciences, Inc. All other service marks, trademarks and tradenames appearing herein are the property of their respective owners.

For Research Use Only. Not intended for use in diagnostic procedures.

R8-50000 Rev. C, Effective Date: 03/17/2025

Cytek Biosciences Inc. | 47215 Lakeview Boulevard, Fremont, CA 94538

Phone: +1 (510) 657-0102 | Fax: +1 (510) 657-0151 | cytekbio@cytekbio.com | www.cytekbio.com